

AllerSnap® Protein Residue Test — Results Record

PEG feeding equipment — post-clean verification swabs Chestnut Tree House Hospice

Month : _____

Area / Kitchen: _____

Tested by: _____

Colour key (30 min incubation at 37 °C):



Pass — green / no colour change



Caution — grey to light purple



Fail — purple

Equipment swabbed	Date	Time	Pass	Caution	Fail	Corrective action taken	Initials
Blender jug & lid							
Blender blade & cap							
Scraper arm & hub							
Blender base / motor unit							
Chopping board (PEG only)							
Allocated PEG preparation knives							
Allocated small equipment							
Storage container for equipment							
Heatproof jug							
Scales							
Sieves & sieve scrapers							

Method: Swab a 10 x 10 cm area of the cleaned, dry surface, snap and squeeze the reagent, mix, then incubate 30 minutes at 37 °C before reading the colour.

Action on Caution: Re-clean the item with Aseptopol E75, disinfect with Pro 20 (1 minute contact), rinse, air-dry and re-test before use.

Action on Fail: Remove the item from use immediately, re-clean and re-test until a Pass is recorded. Report to the supervisor and record on this sheet.

HSE Manager verification (sign & date): _____

Test kit lot no.: _____